User Manual

Version 3.0

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Easy RNA Extraction Kit


Product introduction

This kit can isolate total RNA easily and rapidly from animal or plant tissues (non-polysaccharides and non-polyphenols) and cultured cells. The kit is easy to operate, and it takes only 10 minutes to complete the extraction process. The isolated total RNA is intact with little DNA or proteins, thus it can be used for RNA related assays such as RT-PCR, Northern blot analysis, mRNA isolation, and *in vitro* translation reactions.

Kit Contents

Cat. No.	50 reactions (DR0401050)
Buffer UR1	50 ml
Buffer UR2	15 ml
Buffer CR	10 ml
RNA Elution Buffer	05 ml
HiPure RNA Column	50
2 ml Collection Tube	50
1.5 ml Collection Tube	50
Handbook	1

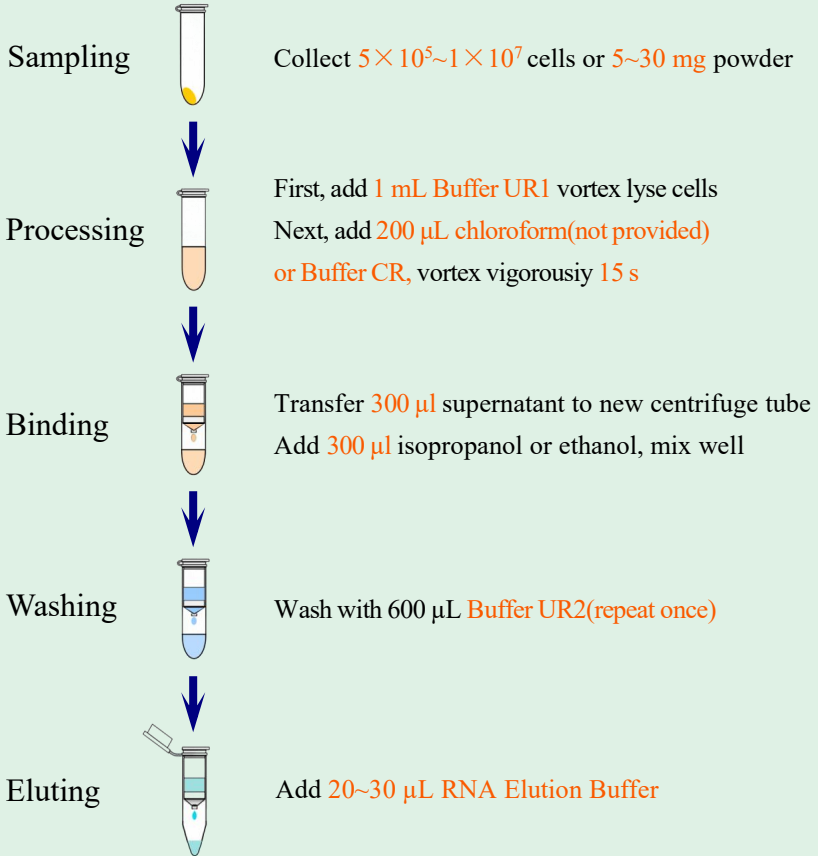Experimental Overview

Treatments of different samples

Cultured cells

Adherent cells: Add 1ml Buffer UR1 onto a 10 cm² petri dish with adherent cells, swirl the reagent around in plate to make sure the dish has been covered with Buffer UR1 (1 ml Buffer UR1 was added in cell culture dish with diameter of 3.5 cm). Repeatedly pipette several times until cells are completely resuspended. Collect the cells with a pipette and transfer them to the 2 ml RNase-free centrifuge tube and the operation proceeds to step 3.

Non-adherent cells: Transfer the cells and media into a centrifuge tube, centrifuge at 8,000 g for 2 minutes at 4 °C, discard the supernatant and avoid disturbing the cell pellet. Add 1 ml Buffer UR1 for every $5 \times 10^5 \sim 1 \times 10^7$ cells. Pipette up and down until cells are fully lysed, and the lysate was transferred to a 2 ml RNase-free centrifuge tube and the operation proceeds to step 3.

Animal and plant tissue sample

The sample was ground to a powder in liquid nitrogen using a mortar and pestle. About 5~30 mg of the tissue powder was weighed in a 2 ml RNase-free centrifuge tube which was precooled with liquid nitrogen, and 1 ml of Buffer UR1 was added, then pipette up and down until cells are fully lysed and the operation proceeds to step 3.

Procedure

1. **Sampling:** Immediately transfer frozen or fresh tissue into mortar, add liquid nitrogen, then crush with pestle to homogenize until powdery. Transfer the tissue powder (about **5~30 mg** powder) into 2 ml RNase-free centrifuge tube.
2. **Lysing:** Add **1 ml Buffer UR1** and then pipette up and down or vortex 15s until cells are fully lysed.
3. **Reacting:** Add **200 μ l chloroform(not provided) or Buffer CR**, vortex and mix 15s, and then centrifuge at maximum speed ($\geq 13,000$ g) for 3-5min.
4. **Binding:** Transfer **300 μ l** top liquid layer to new centrifuge tube without touching middle layer. Add **300 μ l** of **isopropanol or ethanol** for homogenization and mix well. Place a HiPure RNA Column in a provided 2 ml collection tube. To bind RNA, apply **600 μ l mixture** to the HiPure RNA Column, and then centrifuge at 12,000 rpm ($\geq 13,000$ g) for 30 s. Discard flow-through and place HiPure RNA column back into the same collection tube.
5. **Washing:** Add **600 μ l Buffer UR2** to HiPure RNA Column and 12,000 rpm ($\geq 13,000$ g) centrifuge for 30 s. Discard flow-through and place HiPure RNA column back into the same collection tube. Repeat the washing procedure again.

For Research Only

Note: Make sure that 96–100% ethanol has been added to Buffer UR2.

6. **Volatilizing:** Discard the flow-through and centrifuge the HiPure RNA Column for 1 min at 12,000 rpm ($\geq 13,000$ g).
7. **Eluting:** Place HiPure RNA Column into an RNase-free 1.5 ml centrifuge tube. Add **20~30 μ l RNA Elution Buffer***, leave it at room temperature for 1-2 min, centrifuge at 12,000 rpm ($\geq 13,000$ g) for 1 min at room temperature. The liquid in the tube is the solution containing the target RNA.

Note: Heating RNA Elution buffer to 65°C improves elution efficiency.

8. **Testing:** The purified RNA can be checked by agarose gel analysis. The obtained RNA can be used for follow-up experiments immediately or be stored at -80°C.

Storage

The **Buffer UR1 and Buffer CR** should be stored dry at 4 °C while other contents are stored at room temperature, and all caps must be tightly closed to avoid reagent failure. **Add absolute ethanol to Buffer UR2 before use (see bottle label for the detailed volume)**. If the skin or eyes accidentally touch the reagent, especially the **Buffer UR1** which containing organic reagents, rinse it immediately with plenty of water and see a doctor.

Note

All centrifugation steps are performed at room temperature since the addition of Buffer UR1.

Q&A

Q1. Why the extraction rate of RNA is low?

A1. When the RNA extraction rate is low, it can be considered from the following aspects:

- ① The initial sample amount is too little to effectively isolate the RNA, please increase the sample size.
- ② If the initial sample volume is too large, the Buffer UR1 might be not sufficient. Please operate the assay according to the recommended ratio of the kit.
- ③ The polysaccharides and non-polyphenols which present in some plant tissues would form insoluble precipitation with Buffer UR1, causing low extraction rate of RNA.

Q2. Why does the extracted RNA degraded?

A2. When the extracted RNA is degraded, it can be considered from the following aspects:

- ① Tissue used for RNA extraction was not fresh. It is strongly recommended that the samples should be fresh or flash frozen with liquid nitrogen then stored at -80°C .
- ② Too large amount of tissue was used. Add the suggested ratio of Buffer UR1 to the samples.

Q&A

- Q3.** What should I do if the sample size exceeds the amount specified in the manual?
- A3.** This kit is mainly used for small-scale preparation of RNA. When the sample size exceeds the range of the specification, it can be operated through increasing the reagent dosage according to the recommended ratio or dividing into multiple tubes. Too large initial amount will affect the lysis efficiency, resulting in low yield or even experimental failure.

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Zhejiang Easy-Do Biotech CO.,LTD